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comply with the relevant harmonization regulations of the Union as well as other applied directives and standards. The following harmonized standards were used in development and production:

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The declaration of conformity only relates to the design and manufacture of the devices described above in accordance with EU Directive 2014/34 / EU in accordance with the harmonized standards EN ISO 80079-36: 2016 and EN ISO 80079-37: 2016 "constructive safety".

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X = Special conditions for use:

The manufacturer defines the ambient temperature range as $-20^{\circ}\text{C} \leq T_a \leq 60^{\circ}\text{C}$. The intended use of the agitator as well as the explosion protection measures described in the operating instructions in accordance with RL 2014/34 / EU must be observed by the operator. An inductive sensor with ATEX approval for group II, category 2GD for temperature class T6 can optionally be used for speed evaluation.

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TUV®

FULL-SERVICE PROVIDER
FOR DRIVE AND
STIRRING TECHNOLOGY.

TABLE OF CONTENT

| | |
|---|----|
| 1. Preface..... | 26 |
| 2. Safety and responsibility..... | 27 |
| 3. Manufacturer and customer service..... | 28 |
| 4. Transport and storage..... | 28 |
| 5. Intended use..... | 29 |
| 6. Assembly and disassembly..... | 29 |
| 7. Connection overview..... | 32 |
| 8. Commissioning..... | 33 |
| 9. Application..... | 34 |
| 10. Technical data..... | 36 |
| 11. Speed..... | 36 |
| 12. Maintenance..... | 36 |
| 13. Explosion protection..... | 37 |
| 14. Trouble shooting..... | 38 |
| 15. Disposal..... | 39 |
| 16. Certificates..... | 40 |

1. Preface

Thank you for buying a product of PMT mechatronics GmbH. Our products guarantee the highest quality and are manufactured with the latest manufacturing technologies.

In case of any questions, wishes or queries, please do not hesitate to contact us. We will be pleased to assist you.

The operator as well as the user have to familiarize themselves with the device and understand its applications. The operating manual ensures the secure and efficient handling of the device. As the manual is part of the device, it must be kept close to it and readily accessible for personnel.

Additionally, all general safety regulations and the local occupational safety regulations for the application of the device must be observed.

The illustrations in this manual may differ from your device due to the wide variety of construction sizes and equipment packages existing.

All products of PTM mechatronics GmbH are developed and manufactured in accordance with the latest European explosion protection guideline RL2014/34/EU, audited and confirmed by official certification authorities.

2. Safety and responsibility

The user must read this manual before putting the device into operation and adhere to all instructions during operation.

- Repairs must only be conducted by the manufacturer.
- RL 1999/92/EG must be observed (Health protection and safety of employees).
- Please check delivery against your order and delivery note for authenticity and completeness.

All accident prevention regulations and any additional safety regulation of the country, in which the device is operated, must be observed.

| | |
|---|---|
| <p>Compressed air</p>  | <p>Please only use filtered, unlubricated compressed air with 2 to 6 bar and a filtering degree of $\leq 5\mu$.</p> |
| <p>Exhaust air</p>  | <p>The exhaust air can be discharged in a targeted manner via the supplied exhaust air hose.</p> <p>If this is not desired, only the exhaust air silencers optionally available from PTM may be used.</p> |

3. Manufacturer and customer service

PTM mechatronics GmbH
Gewerbepark 1
82281 Egenhofen
Germany

Phone: +49 8134 - 25797 - 0
Fax: +49 8134 - 25797 - 99
Email: info@ptm-mechatronics.com
Website: www.ptm-mechatronics.com

Please contact our customer service before shipping returns or repair devices. Please indicate the device number.

4. Transport and storage

The device is delivered in a carton package which protects it against any damages. Please make sure that the package is not damaged.

For storage, please observe the following instructions:

- Store dry
- Avoid shocks
- Storage temperature: 10 to 40 degrees Celsius
- Do not storage in damaged packaging

Please inspect all parts and packaging regularly in case of long storage. Please replace damaged packaging if necessary. After unpacking the device, please check for completeness and transport damages.

In case of any visible transport damages, please follow the following instructions:

- Do not accept delivery or only under reserve.
- Note down extent of damage on transport documentation or delivery note of carrier.
- Initiate complaints processing.

5. Intended use

Agitators and stirrer drives are used for the drive of stirrer elements, agitator shafts and axes, for example for the stirring and mixing of liquids and powders in open or closed, pressureless containers. Both, agitators and drives, are referred to as „agitators“ in this manual. For a safe operation, agitators must be mounted firmly to a housing, to a tripod, to the cap of a container or a lifting station. They are powered by compressed air. Agitators are available in anodised aluminium or stainless steel (1.4404). Agitator shafts and elements are made of stainless steel (1.4404).

Agitators are available in different versions. For more information, please refer to: www.ptm-mechatronics.com

6. Assembly and disassembly

6.1. General safety instructions

- Impacts involving rust and light metals with their alloys may cause ignitable sparks.
- The use of tools with corroded surfaces is prohibited.
- When choosing materials and fixing accessories, please consider corrosion, abrasion and other interactions.
- Ensure potential equalisation.
- Consider tightening torques correspondingly.
- Ensure correct installing of stirrer elements.

All accident prevention regulations and possible additional safety regulations of the respective country must be observed.


6.2. Mounting of impellers

The agitator shafts (round or hexagonal) are attached to the laboratory agitator using the quick-release chuck.

When mounting impellers, please always observe a correct alignment as described as follows.


6.2.1. eco-Disc

The larger diameter of the eco-Disc cups must point in the direction of rotation.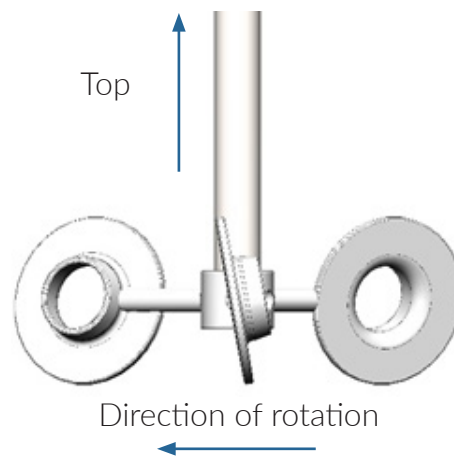


6.2.2. eco-Cup

The larger diameter of the eco-Cup cups must point in the direction of rotation.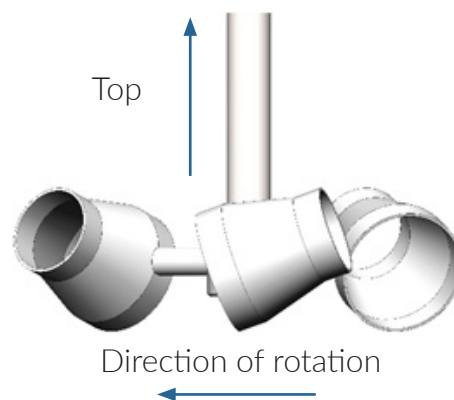


6.2.3. Dissolver disc

For the correct alignment, please follow the mark on the dissolver disc.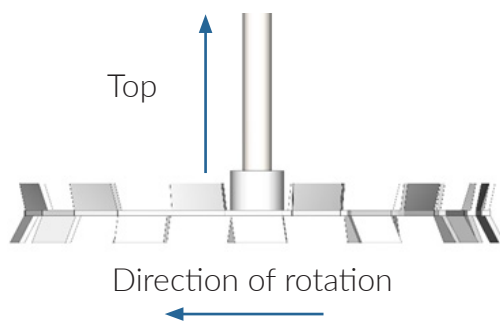


6.2.4. Helical ribbon impeller

Before mounting, please make sure that the impeller is not damaged or bent out of shape.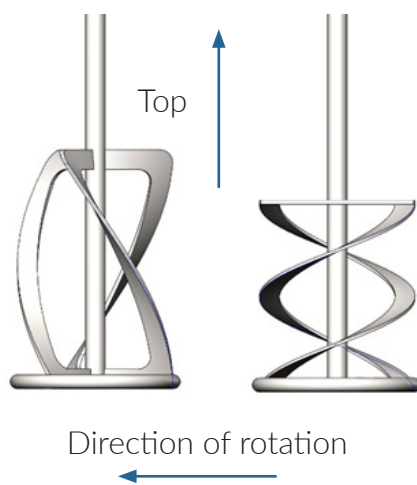

7. Connection overview

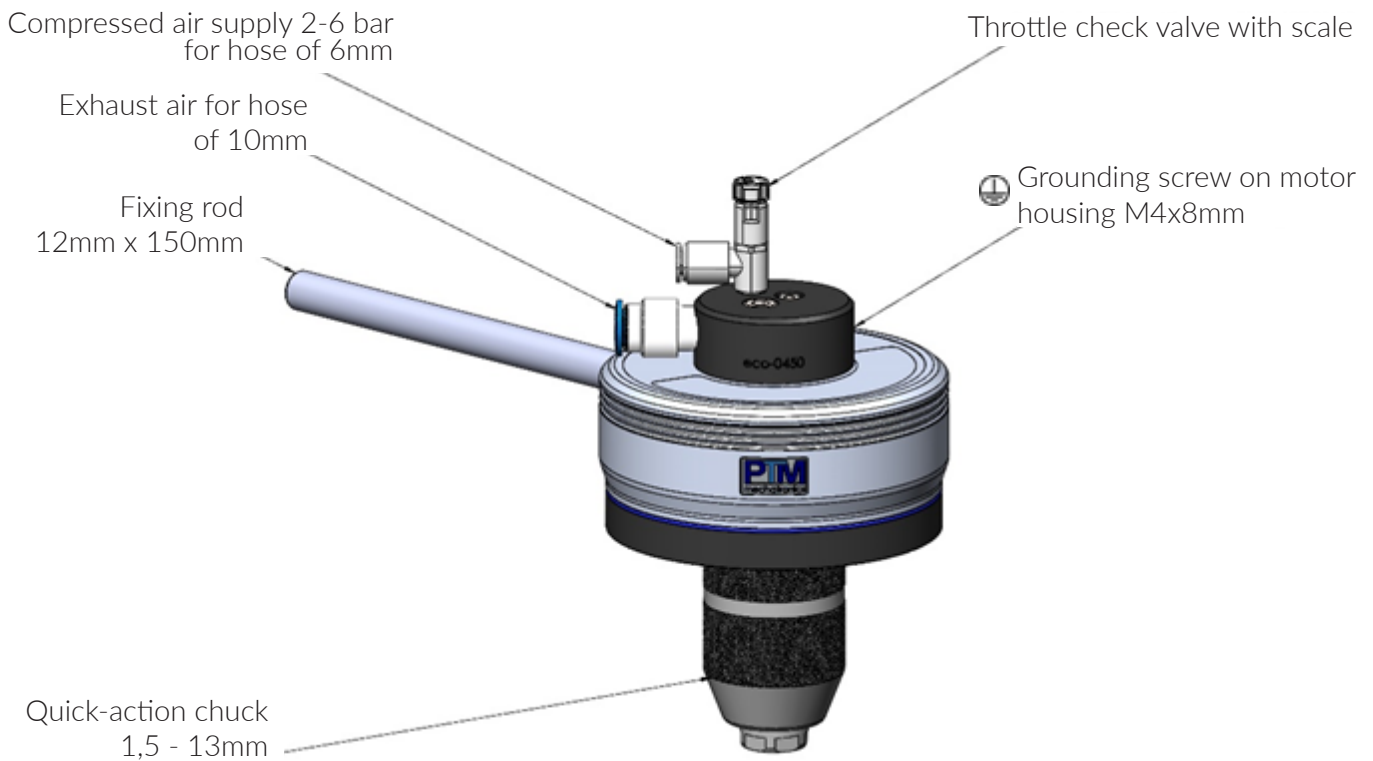Fixing of shaft with quick-action chuck:

- Open chuck by gripping retaining ring and turning it to the left
- Push agitator shaft into open chuck
- Close chuck by gripping retaining ring and turning it to the right
- Chuck snaps when closed to secure impeller adequately

Handling:

- Turn off throttle check valve
- Connect grounding ⚡ to motor housing
- Connect air supply
- Connect exhaust hose and fix it with a downward gradient
- Attach agitator shaft with quick-action chuck and impeller
- Mount agitator with fixing rod, for example on tripod
- Only use agitator in filled containers
- Set speed by means of throttle check valve

Grounding can take place via the grounding screw or via the air connection and an electrically conductive air supply hose. This supply hose must be grounded.

8. Commissioning

8.1. General

To guarantee the proper function of the agitators and ensure the longest possible service life, the following instructions must be observed when putting the devices into operation:

- All packaging remains must be removed thoroughly.
- All connections must be fixed accurately.
- For supply air, please make sure to use the hose diameter required according to supply air throttle valve.
- All main air lines must have a gentle gradient downward of at least 1% in the direction of flow, thus enabling draining of condensation water at the lowest point by means of a valve (manually or automatically).
- In case of horizontal main pipes, always connect junctions on the upper side, in case of vertical main pipes, do not connect junctions at the lowest point, to prevent condensation water from flowing into the junction.
- Use maintenance units with an air filter with a maximum filtering degree of 5 μ Counter nut
- Install maintenance unit with an air connection of at least G 1/4" in close proximity to the agitator.
- The air filter must be maintained regularly. Please clean the filter and let the collected condensation water drain off.
- All agitators provide their best performance at 6 bar (85psi). Should the pressure in the compressed air net exceed 6 bar, a pressure regulator must be used.
- Maximum speed must not be exceeded.
- Use the exhaust hose delivered with your device.
- Connect exhaust hose with a downward gradient.
- Do not close the exhaust hose. Do not use silencers.

All agitators comprise a housing of protection class IP 67, which guarantees the necessary IP protection. Nevertheless, the date for cleaning the device to remove dust must be determined by the operator.

8.2. Verifying of explosion protection

Before commissioning in hazardous environments, please verify:

- Is the agitator explosion-protected?

- Is the unit explosion-protected?
- Is the equipotential bonding provided?
- Are conductive hoses and connections used?

8.3. Compressed air

Function and service life of agitators largely depend on quality and pressure of the compressed air used. High humidity or dirt particles destroy the agitator. A high percentage of condensate leads to rust inside of the agitator. This is why it is important to observe the following:

- Agitators are only admitted for the operation with compressed air of not less than quality class 3 according to DIN ISO 8573-1 which was generated and processed outside of the Ex area.
- The use of liquids or gases to operate the devices does not comply with the intended use.

9. Application

9.1. General

- Agitators are only admitted for the appropriate and intended use. In the event of violation, warranty and manufacturer responsibility expire instantly.
- Manipulation of all kinds leads to expiry of conformity, warranty and manufacturer responsibility.
- Agitators may only be operated with a maximum pressure of 6 bar.
- Speed ranges must be adhered to (please refer to section 27).
- The environmental conditions specified in this manual must be observed.
- The operator must ensure the correct mounting of agitator shaft and stirrer element.
- The operator must ensure that no foreign objects get into the container which may cause sparks during operation.

9.2. Application in Ex area

The agitators can be used in the following hazardous areas:

- in zone 1
- in zone 21
- in zone 2
- in zone 22

Please observe the following instructions for the application in Ex areas:

- The agitator may be put into operation in categories 2 and 3 (zone 1 or 21, zone 2 or 22) only by specialists.
- Exhaust air is allowed to escape in ATEX zone 1 and 2.
- All closures must be attached again after opening and closing.
- Openings for supply and exhaust air must not be closed.
- In case of discharge of electrostatically charged parts, ignitable sparks may be caused. This is why hose and bundles of hoses should not exceed a maximum outer diameter of \varnothing 20 mm.
- For operation, an adequate equipotential bonding must be established with the grounding screw on the housing.
- Agitators must be protected against objects falling down.
- Operation is only permitted with closed and intact housings.
- Operation of damaged devices is forbidden.
- Number and variety of removable connections must be limited to a minimum. Hose connections must be held as short as possible. Mechanical tensions must be avoided.
- Protection class IP 67 is sufficient for housing and side parts. An additional protection against humidity is not necessary.
- In hazardous areas, only accessories may be used that fulfil all requirements of the latest European guidelines and the respective national law.
- Direct dangerous electrostatic discharges on the agitators are forbidden. Discharges like that normally cannot be caused by humans but require a compressed air nozzle or the like.
- Operation in dust Ex areas is permitted in case of non-conductive dusts with an ignition energy of 3 mJ or higher.
- The operator must ensure adequate lighting protection measures.
- Stirring in non-conductive containers requires a conductivity of the liquid of $> 10000\text{pS/m}$. Otherwise, the operator must inert the container or add conductivity additives.
- Pass-through operations are permitted provided that all guidelines of TRGS 727 are observed. Stirring container and stirrer element must be conductive and grounded. The liquid stirred must have a high conductivity. Agitator speed must be reduced to avoid squirting or spraying of the liquid. These guidelines must be fulfilled by the operator.
- To avoid contact between shaft and container wall, attention must be paid to a sufficient spacing between shaft and container wall when dimensioning the shaft. Shafts and stirrer elements that are purchased from PTM mechatronics always fulfil these requirements. Otherwise, the operator must ensure an adequate dimensioning.
- The operator must prevent any development of heat due to chemical reactions between two substances.

10. Technical data

| | |
|--|--|
| Connection compressed air | 2.0 bar – 6.0 bar |
| Ex marking gas | II 2 G Ex h IIC T5 Gb X |
| Ex marking dust | II 2 D Ex h IIIC T100°C Db X |
| CE marking | CE |
| Liquid temperature | max. +50° Celsius |
| Ambient temperature | -10° to + 60° Celsius |
| Speed (at 6 bar, 150 ¹ /min) | 50W to 230W (depending on construction size) |
| Air consumption (at maximum load) (construction size) | 25 l/min to 500 l/min (depending on construction size) |

11. Speed

| | |
|---------------------------------------|---|
| Stirrer drives without transmission: | 0 ¹ /min -300 ¹ /min |
| Stirrer drives with 1:2 transmission: | 0 ¹ /min - 600 ¹ /min |

Speed ranges specified above are subject to a tolerance of +/- 10%.

12. Maintenance

12.1. General

When maintaining agitators that are operated in hazardous areas, the operator must ensure that all technical regulations for operational safety, the Industrial Safety Regulation and all requirements of the respective country are adhered to.

- Maintenance measures must only be conducted by PTM mechatronics GmbH or experts especially trained by PTM mechatronics GmbH.
- Components must only be replaced by original spare parts that are certified for the application in Ex areas. The same applies for lubricants and auxiliary materials.
- Devices must be maintained and cleaned regularly. Intervals are to be determined by the operator considering the environmental circumstances on site, for example dust deposits of approximately 0.5 - 1 mm.
- After maintenance, all removed barriers and warnings must be reattached to their former position.
- Dust deposits in gaps of movable parts can potentially heat up. Therefore, it is necessary to clean those gaps regularly.
- Agitators must only be cleaned outside of hazardous areas.
- Wear parts include pistons, ball bearings, seals and shaft seals.

12.2. Pistons

The pistons of radial piston motors are wear parts. For a long service life it is very important to examine the motor and its pistons regularly. We recommend an immediate examination of the motor and its pistons in case the motor sound changes or increases or the motor loses performance.

We recommend the following examination intervals:

| | |
|-----------------------------------|-----------------|
| Year-round, three-shift operation | every 6 months |
| Year-round, two-shift operation | every 12 months |
| Year-round, one-shift operation | every 24 months |
| Irregular, left/right operation | every 48 months |

Pistons should be replaced immediately in case of wear to avoid major damages.

12.3. Bearings

Bearings must be replaced after a service life of 35.000 hours at most when the agitator is used in hazardous areas. To analyse speed an inductive sensor of ATEX category 2 can be used optionally.

12.4. Seals and shaft seals

Seals and shaft seals are wearing parts. The factors that have an unfavorable influence on wear are the operating hours and the speed. Other factors that promote wear include environmental conditions such as temperature, dust, solvents and vapors.

The operator is responsible for assessing wear. If necessary, the seals and shaft seals must be replaced.

13. Explosion protection

13.1. Basics

Due to its planned application in hazardous areas, the following EU guideline is decisive for the development of the agitator:

European explosion protection guideline: 2014/34/EU

13.2. European guidelines, laws and provisions

2014/34/EU: European explosion protection guideline

13.3. Marking according to guideline 2014/34/EU and normative requirements

Marking of non-electric equipment is:

 II 2 G Ex h IIC T5 Gb X
 II 2 D Ex h IIIC T100°C Db X

13.4. Explosion protection provisions

For devices operated in hazardous areas, the operator must observe a number of provisions. The following list gives an overview of the most important provisions.

Within the European Union the following provisions apply:

- Directive 1999/92/EC regarding minimum regulations for improving health safety and safety of employees who can be endangered by hazardous atmospheres.
- DIN EN 60079-0 regarding electrical equipment for hazardous areas – General Provisions
- DIN EN 60079-10 regarding hazardous areas - Section 10-1: Classification of hazardous areas
- DIN EN 60079-14 regarding electrical equipment for potentially gas explosive areas - Section 14: Electrical systems in hazardous areas
- DIN EN 1127-1 regarding explosive atmospheres – explosion protection - Section 1: Principles and methods
- DIN EN ISO 80079-36 Non-electrical devices for application in hazardous areas – Section 1: Principles and requirements
- DIN EN ISO 80079-37 Non-electrical devices for application in hazardous areas – Section 5: Protection through constructive safety „c“
- TRGS 727 Technical rules for hazardous substances „Prevention of danger of ignition due to electrostatic charge“

Additionally, further national and international regulations and guidelines may apply.

13.5. Zone classification for hazardous areas

Hazardous areas are areas in which hazardous atmospheres may occur to a dangerous extent due to local and operational circumstances. They are subdivided in zones for combustible gases, vapours or mist:

- Zone 0/20 – includes areas in which hazardous atmospheres are present constantly or over the long term.
- Zone 1/21 – includes areas in which hazardous atmospheres can be expected to occur occasionally.
- Zone 2/22 – includes areas in which hazardous atmospheres can be expected to occur only seldom and even then, only for a brief period.

14. Trouble shooting

In case the agitator does not perform adequately, please check the following points:

- Is the air pressure sufficient? (2-6 bar / 85 psi)?
- Is the air pressure stable?

- Does the air system provide a sufficient amount of air?
- Is the air filter of the maintenance unit clean?
- Are there any leaks between maintenance unit and motor?
- Are there any cross-section constrictions in the exhaust line?

Should none of the above mentioned defects match, the agitator must be sent in to the manufacturer for examination.

Devices that are operated in hazardous areas must not be modified. Repairs must only be conducted by the manufacturer or by experts especially trained by the manufacturer.

15. Disposal

Disposal of packaging and used parts must comply with the regulations of the respective country in which the device is operated.

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Zertifikatsinhaber: PTM mechatronics GmbH
TÜV SÜD Product Service GmbH • Zertifizierstelle • Ridlerstraße 65 • 80339 München • Deutschland