Schnellspannsystem „eco-Quick V2“

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Quick-Release System „eco-Quick V2“

eco-Quick V2 is a quick-release system designed to make changing agitator shafts easier and faster.

Always ensure that the agitator drive is secured against being switched on again before replacing the agitator shaft.

Before connecting the agitator shaft with the eco-Quick V2, make sure that both the agitator shaft and the eco-Quick V2 have been thoroughly cleaned and are free from damage.

Figure 1 shows the sequence of steps for installing the agitator shaft with eco-Quick V2:

6. With one hand, pull down the clamping sleeve and the splash guard sleeve of the eco-Quick V2 and then hold them in the lower position (steps '1' and '2').
7. With your free hand, insert the cleaned agitator shaft with its upper end into the eco-Quick V2 (steps '3' and '4').
8. Twist the agitator shaft against the clamping sleeve and the splash guard sleeve (while keeping both sleeves held firmly) until the clamping sleeve and the splash guard sleeve automatically snap back upward (step '4').
9. The agitator shaft is now secured against falling out (step '5').
10. To remove the agitator shaft, perform steps '1' and '2', while holding the agitator shaft firmly with your free hand.

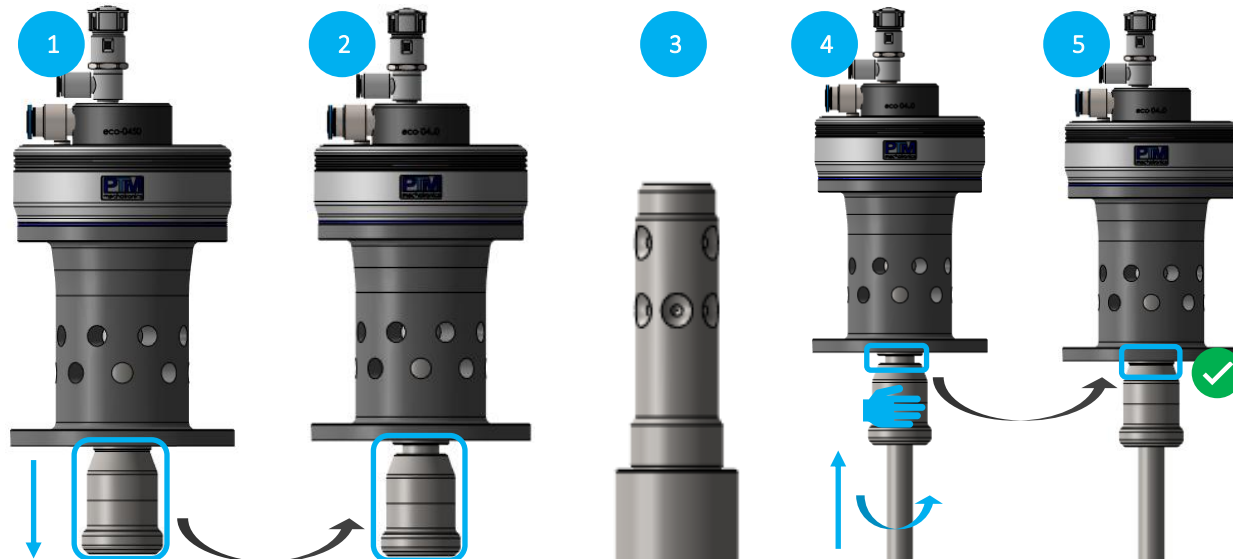
Figure 1: Assembly of the agitator shaft using eco-Quick V2

Quick-Release Systems „eco-Quick V3“ & „eco-Quick V4“

eco-Quick V3 and eco-Quick V4 are quick-release systems designed to make changing agitator shafts easier and faster.

Always ensure that the agitator drive is secured against being switched on again before replacing the agitator shaft.

Before connecting the agitator shaft with the eco-Quick V3 or eco-Quick V4, make sure that both the agitator shaft and the eco-Quick V3 or eco-Quick V4 have been thoroughly cleaned and are free from damage.

Figure 2 shows the sequence of steps for installing the agitator shaft with the eco-Quick V3 (installation of the agitator shaft with the eco-Quick V4 is carried out in the same way as with the eco-Quick V3).

6. With one hand, pull down the splash guard sleeve of the eco-Quick V3/V4 and hold it in the lower position (steps '1' and '2').
7. With your free hand, insert the cleaned agitator shaft with its upper end into the eco-Quick V3/V4 (steps '3' and '4').
8. Twist the agitator shaft against the sleeve of the eco-Quick V3/V4 (while holding the sleeve) until the splash guard sleeve automatically snaps back upward (step '4').
9. The agitator shaft is now secured against falling out (step '5').
10. To remove the agitator shaft, perform steps '1' and '2', while holding the agitator shaft with your free hand.

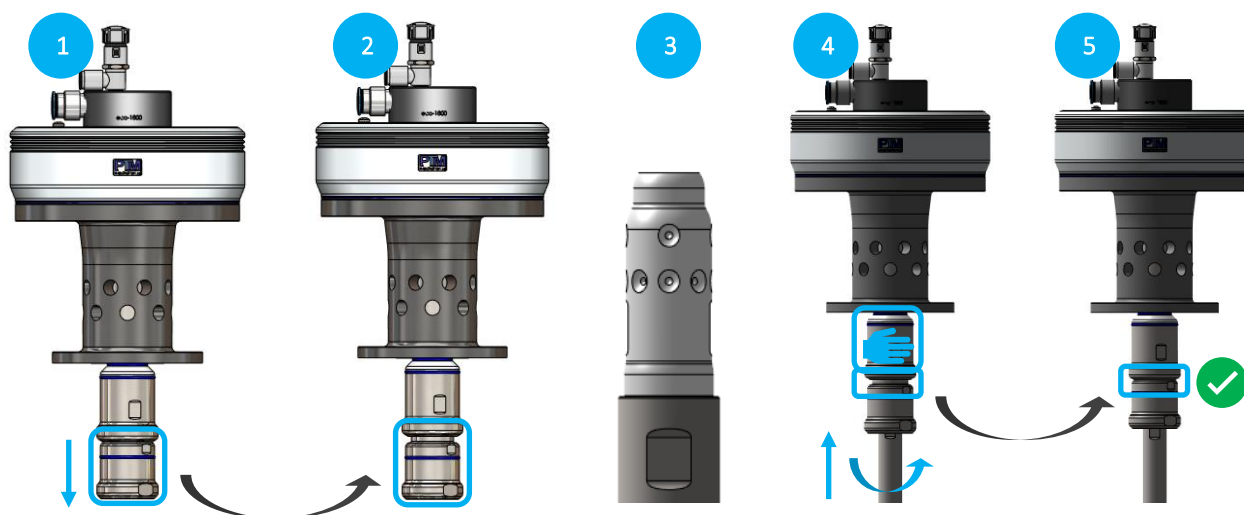
Figure 2: Assembly of the agitator shaft using eco-Quick V3 or eco-Quick V4

Version des Dokuments / Document version: V1